



VeraCode® ADME Software v1.0.1.4 Release Notes

This release note describes new features and changes since the initial VeraCode ADME Software commercial release.

1. Version 1.0.1.4

Note: Uninstall VeraCode ADME Software v1.0.1.3 before installing the VeraCode ADME Core Panel Kit v1.1.8.6 and software v1.0.1.4.

- 1.1. Updated the special case default wild-type star call for NAT1 and NAT2 from *1A/*1A to *4/*4, because the scientific reference changed.

2. Version 1.0.1.3

- 2.1. Changed the STC scoring algorithm to be more tolerant if wells have a zero bead population for STC assays.
- 2.2. Excluded assays with zero beads from the input to normalization algorithm.

3. Version 0.9.4.0 (Initial Commercial Release)

FOR RESEARCH USE ONLY

© 2012–2013 Illumina, Inc. All rights reserved.

Illumina, *illumina*Dx, BaseSpace, BeadArray, BeadXpress, cBot, CPro, DASL, DesignStudio, Eco, GAllx, Genetic Energy, Genome Analyzer, GenomeStudio, GoldenGate, HiScan, HiSeq, Infinium, iSelect, MiSeq, Nextera, NuPCR, SeqMonitor, Solexa, TruSeq, TruSight, VeraCode, the pumpkin orange color, and the Genetic Energy streaming bases design are trademarks or registered trademarks of Illumina, Inc. All other brands and names contained herein are the property of their respective owners.