Local Run Manager v4 への TruSeq DNA CD Index Kit の追加手順

<事前準備>

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The image shows the 'Local Run Manager' web interface. The top navigation bar includes 'Local Run Manager', 'ISEQ', 'ISEQ100', 'RUN DASHBOARD', and 'TOOLS'. The user is logged in as 'Admin User'. The main content area is titled 'Index & Library Prep Kits' and features a 'LIBRARY PREP KIT' tab selected and highlighted with a red box. There are buttons for 'Download Template' and 'Add Index Kit'. Below the tab is a table with the following data:

NAME	DESCRIPTION	SUPPORTED LIBRARY PREP KITS	VERSION
TruSeq Custom Amplicon	TruSeq Custom Amplicon Index Kit	TruSeq Custom Amplicon	1
TruSeq Custom Amplicon Low Input	TruSeq Custom Amplicon Low Input Index Kit	TruSeq Custom Amplicon Low Input	1
TruSeq DNA CD Indexes 96 Indexes	TruSeq DNA CD Indexes 96 Indexes Index Kit	TruSeq DNA Exome Enrichment, TruSeq DNA PCR-Free, TruSeq Nano DNA	1
TruSeq DNA Methylation	TruSeq DNA Methylation Index Kit	TruSeq DNA Methylation	1
TruSeq DNA Single Indexes Set A	TruSeq DNA Single Indexes Set A Index Kit	TruSeq DNA Exome Enrichment, TruSeq DNA PCR-Free, TruSeq Nano DNA	1
TruSeq DNA Single Indexes Set A B	TruSeq DNA Single Indexes Set A B Index Kit	TruSeq DNA Exome Enrichment, TruSeq DNA PCR-Free, TruSeq Nano DNA	1
TruSeq DNA Single Indexes Set B	TruSeq DNA Single Indexes Set B Index Kit	TruSeq DNA Exome Enrichment, TruSeq DNA PCR-Free, TruSeq Nano DNA	1
TruSeq Exome - 8 rxn x 3 6 9 plex	TruSeq Exome - 8 rxn x 3 6 9 plex	TruSeq DNA Exome Enrichment	1
TruSeq Exome - 8rxn x 12plex	TruSeq Exome - 8rxn x 12plex Index Kit	TruSeq DNA Exome Enrichment	1
TruSeq Methyl Capture EPIC	TruSeq Methyl Capture EPIC Index Kit	TruSeq Methyl Capture EPIC	1

At the bottom of the table, there is a pagination control showing '7' selected and 'Showing 81-70 of 83 item(s)'.