

Install Notes

cBot v2.0.29

Recipe Installer v2.0.5

For cBot

Introduction

These instructions detail how to install cBot software version 2.0.29. These update instructions apply when upgrading from any version of cBot software v1.1 or higher.

NOTE: installing the cBot 2.0.29 software as described below will also update and install all recipes.

To confirm the current version of cBot software, select "About" from the Menu tab and the current version will be displayed:

For full information on updated features in cBot 2.0.29 software, please refer to the cBot 2.0.29 Release Notes prior to installation.


I. Upgrading from pre-cBot 1.3 to cBot 2.0.29

For instructions on how to upgrade your cBot software if you are on a version earlier than 1.3, please contact your FAS/FSE or Illumina Product Support.


II. Upgrading cBot 1.3 or later to cBot 2.0.29

Using software v1.3 or later, you can upgrade instrument software from the configuration screen using a USB flash drive. To download the cBot 2.0.29 software, go to http://support.illumina.com/sequencing/sequencing_instruments/cbot/downloads.ilmn and scroll down to cBot 2.0.29 Software. Download the 2.0.29 package and copy it onto a USB flash drive. Do not unzip the package.


1. Insert the USB flash drive containing the new software version installer into one of the two USB ports on the front of the instrument.
NOTE: The installer must reside in the root directory of the USB flash drive, not in a folder. Ensure only the most recent version of installer is present.
2. Touch Menu in the upper-left corner of the screen, and select Configure. The keyboard appears.
3. Log in using the default password **admin** and touch "Enter" to close the keyboard. The configure screen opens as shown below:

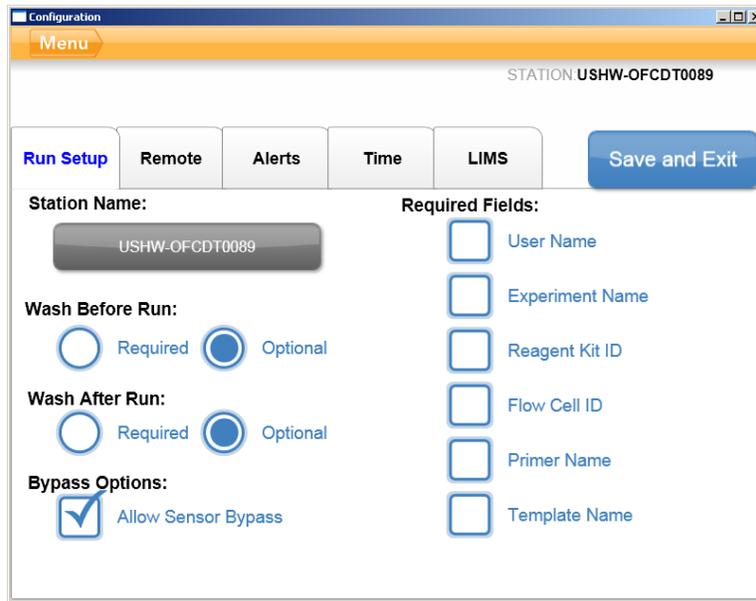

4. Touch Menu in the upper-left corner of the screen, and select Upgrade:

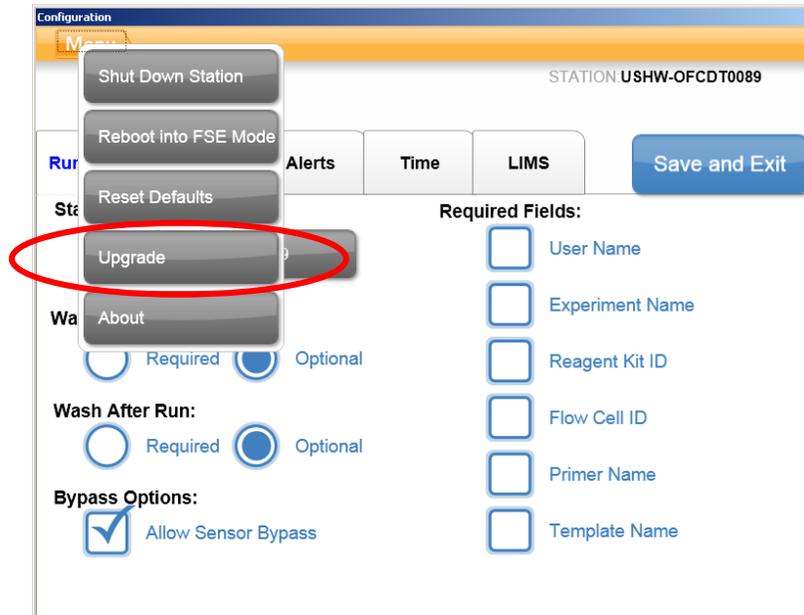

5. If a valid cBot installer cannot be located on either of the 2 USB drives, the message below appears. Either insert a valid USB key containing the cBot upgrade installer and click OK to try again, or cancel out of the upgrade operation.

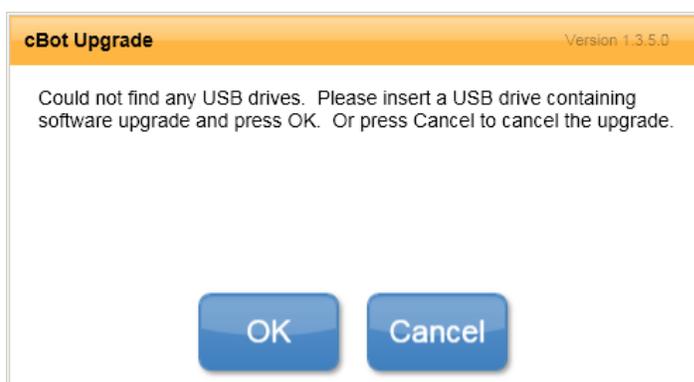

6. Assuming a valid cBot installer is located on either USB drive, a dialog box appears with one of the following messages:

Message	Action
The software installer version is greater than the version currently installed on the cBot	Touch OK to proceed with the installation of the newer version.
cBot cannot find a valid software installer	You can either insert a valid cBot upgrade and touch OK to try again, or Cancel to abort the upgrade.
The software installer version is equal or lower than the version currently installed on the cBot	Touch Cancel to abort the upgrade, or OK to proceed with installation of a previous version.

WARNING: The USB flash drive must remain in the USB slot for the duration of the upgrade process until the instrument reboots automatically. Do not interact with the system during the upgrade.

Upon completion of the upgrade, the cBot instrument reboots automatically with the new software version. The reboot process takes about ten minutes to complete.

7. When the cBot reboot is complete and the login screen appears, you can remove the USB flash drive.
8. Following installation, please confirm in the About box (as detailed in the introduction) that the software has been updated to version 2.0.29. Note that the cBot 2.0.29 installation will update with the recipes required for use with HiSeq X v2 HD cluster and HiSeq X v2 rehyb workflows.

III. Upgrading recipes in cBot 2.0.29

NOTE: installing the cBot 2.0.29 software as described above will also update and install all recipes, so completing the section below will not be necessary in most cases. Upgrading to recipe installer v2.0.5 without upgrading to cBot 2.0.29 software will not support HiSeq X v2 barcode masks.

The recipe installer and these instructions are provided in case there is a need to update recipes independent of updating cBot software. Using the cBot software v2.0.29 package, you can upgrade recipe versions independent of software upgrades using a USB flash drive containing the recipe installer.

The cBot 2.0.29 software update will automatically hide or archive v7 GA recipes. If these are still needed, move the v7 recipes from C:\Illumina\cBot\bin\Recipes\Hidden to C:\Illumina\cBot\bin\Recipes.


1. Insert the USB flash drive containing the Recipe Installer 2.0.3 .msi into one of the two USB ports on the front of the instrument.
NOTE: The installer must reside in the root directory of the USB flash drive, not in a folder. Ensure only the most recent version of installer is present.
2. Touch Menu in the upper-left corner of the screen, and select Configure
3. Log in using the default password **admin** and touch enter to close the keyboard. The configure screen opens as shown below:

The screenshot shows the 'Configuration' window for station 'USHW-OFCDT0089'. It features a 'Menu' button in the top left and a 'Save and Exit' button in the top right. The main area is divided into several sections:

- Station Name:** A text field containing 'USHW-OFCDT0089'.
- Wash Before Run:** Two radio buttons, 'Required' (unselected) and 'Optional' (selected).
- Wash After Run:** Two radio buttons, 'Required' (unselected) and 'Optional' (selected).
- Bypass Options:** A checked checkbox for 'Allow Sensor Bypass'.
- Required Fields:** A list of six checkboxes, all unselected: 'User Name', 'Experiment Name', 'Reagent Kit ID', 'Flow Cell ID', 'Primer Name', and 'Template Name'.

Navigation tabs at the top include 'Run Setup' (active), 'Remote', 'Alerts', 'Time', and 'LIMS'.

4. Touch Menu in the upper-left corner, and select Upgrade Recipes:

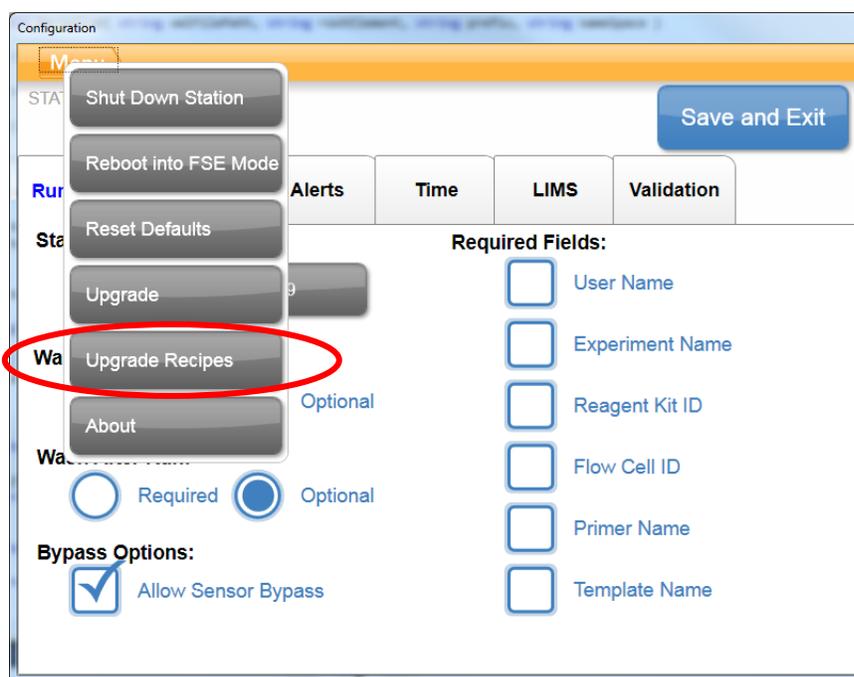

WARNING: The USB flash drive must remain in the USB slot for the duration of the upgrade process until the instrument reboots automatically. Do not interact with the system during the upgrade.

Upon completion of the upgrade, the cBot reboots automatically with the new recipes. The reboot process takes about ten minutes to complete.

5. When the cBot reboot is complete and the login screen appears, you can remove the USB flash drive.

IV. Known issues

On a small subset of instruments it has been observed that the installer appears to run and complete but the software does not get updated and remains on the 1.4 version.

If this occurs please download the old cBot1.4 software and performing a repair install of that software first, followed by rerunning the new cBot 2.0.29 software installer, as shown in the instructions above.

Installation Troubleshooting:

Please contact Illumina Technical Support if any other issues arise during installation.