



# **iSeq™ 100 System Release Notes**

**Suite Installer v1.3.0**

**iSeq™ 100 Control Software Server v1.3.0**

**iSeq™ 100 Control Software Client v1.3.0**

**Firmware (FW) v1.3.0**

**Real-Time Analysis (RTA) v2.12.5**

**Universal Copy Service (UCS) v1.5.4**

**Local Run Manager v2.1.2**

**17-AUG-2018**

## Introduction

These release notes detail new features, improvements, and known issues for the iSeq 100 Sequencing System. For more information on the iSeq 100 System, see the *iSeq 100 Sequencing System Guide* (document # 1000000036024) on the Illumina website.


The software package includes:

- Suite Installer v1.3.0
- iSeq™ 100 Control Software Server v1.3.0
- iSeq™ 100 Control Software Client v1.3.0
- Firmware v1.3.0
- Real-Time Analysis (RTA) v2.12.5
- Universal Copy Service v1.5.4
- Local Run Manager v2.1.2

## I. Suite Installer v1.3.0

### KNOWN ISSUES

- The following prompt will appear during software install.  
Select **Yes**. If No is selected, contact Illumina Technical Support. This prompt enables Software Restriction Policies (SRP). See the *iSeq 100 Sequencing System Guide* for more information.

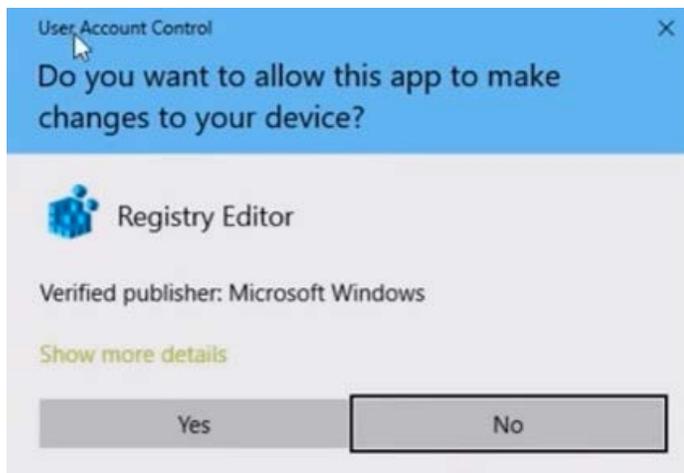

## II. iSeq Control Software Client v1.3.0

### NEW FEATURES

- The "Save thumbnail" option now includes additional images to aid in run debugging.
- Average Q30 for Read 1, Read 2, and projected output is written to run completion file.

- Added ability to access Local Run Manager through the control software drop-down menu
- Enabled editing of Universal Copy Service account via the Network Configuration tab in System Settings.
- Enhancements and robustness added to Pre-Run Check for imaging, storage space, and transfer checks.
- Updates to system check specifications for reliability and robustness.

#### BUG FIXES

- Eliminated intermittent failures to connect to embedded system during instrument initialization.
- BaseSpace™ Sequence Hub communication error is displayed on the correct screen during first time setup.
- Process Management page wording updated upon run deletion failure.
- New runs no longer prevented due to errors from previous run.
- Multiple updates to error messages for UCS connection loss, software update failures, cartridge loading errors, and system check failures. These changes improve clarity and suggested corrective action.
- Saving of system settings supported despite BaseSpace Sequence Hub communication error.
- User interface updates.
- Resolved error generated when clicking Q30 values on run screen.

#### KNOWN ISSUES

- Using a sample sheet on a mapped network drive in manual mode results in a run failure. To avoid this failure, move sample sheets to the local hard drive before starting a run.
- Minimizing control software during a sequencing run might result in pre-run checks or run failure.
- Immediately launching the System Check screen after launching control software might freeze the user interface. Allow 3–5 seconds after initialization before launching system checks.

### III. Universal Copy Service v1.5.4

#### NEW FEATURES

- Enhanced robustness of file copying and upload process

#### KNOWN ISSUES

- No known issues.



## **IV. Local Run Manager v2.1.2**

### **NEW FEATURES**

- Added storage space availability check.
- Run setup unable to proceed when Local Run Manager does not have access to the user designated run output folder.

## **V. Server v1.1.0**

- No known issues or new features.

## **VI. Real-Time Analysis v2.12.5**

- No known issues or new features.

## **VII. Firmware v1.3.0**

- No known issues or new features.