



TruSight Tumor 15 LRM Customer Release Notes
Part Number: 200003521
Ver: 00
Effective Date: 06-Aug-2021 to UP (NONE)
Release Status: Released
Release Date: 01-Jul-2021
IP Classification: --

Page **1** of **2**

TruSight Tumor 15

LRM Module

Release Notes

V2.1.0

For TruSight Tumor 15 Reagents



Introduction

These Release Notes detail the key features and known limitations for the TruSight Tumor 15 v2.1.0 Local Run Manager module.

NEW FEATURES:

- Overhauled the PDF reporting infrastructure, using now Headless Chrome to create PDF reports. This new infrastructure is more stable and fixes a previous bug where some rows would be partially cut out.
- Updated the Illumina Annotation Engine to version 2.0.11
- Added support for LRM version 3.0 (MiSeq) and LRM version 2.4 (MiniSeq).
- LRM version 1.3 (MiniSeq) is no longer supported.
- Bcl2FastQ index mismatch is now set to zero to address an issue with index mis-assignment. A full explanation of this issue can be found in PQN2020-1086.

DEFECT REPAIRS:

- Fixed issue where requeuing an analysis would skip the Edit Setup stage if performed by an admin user.
- Fixed issue where a null value for Consequence resulted in the HTML report in LRM not being populated.
- Updated installer so that installation and/or uninstallation does not fail in later LRM versions when not specifically run from an administrator window.