

# MiniSeq Software Release Notes

**MiniSeq System Suite v2.1.0**

**MiniSeq Control Software v2.1.0**

**MiniSeq Service Software v2.1.0**

**MiniSeq Recipes v2.1.0**

*For MiniSeq Systems*

## Revision History

Revision	Date	Description of Change
00	21/SEP/2016	Initial release
01	17/FEB/2020	MiniSeq v2.0 software changes
02	24/AUG/2020	MiniSeq v2.1 software changes

## Introduction

These Release Notes detail new features, improvements, and issue fixes for the for the Illumina MiniSeq Software. For more information on the MiniSeq Sequencing System and associated software, see the MiniSeq Sequencing System page on the Illumina Support Site.

This release is only compatible on Microsoft Windows 10. If your system is running on Microsoft Windows 7 and would like to upgrade to the latest Microsoft Windows 10, reach out to your local sales representative or support team to schedule an upgrade.

Software Application	Prior Version	New Version
MiniSeq System Suite Installer	2.0.0	2.1.0
MiniSeq Control Software (MNCS)	2.0.0	2.1.0
MiniSeq Service Software (MNSS)	2.0.0	2.1.0
MiniSeq Recipe Installer	2.0.0	2.1.0
MiniSeq Real Time Analysis (RTA)	2.11.4	2.11.4
Universal Copy Service	1.6.3	1.6.3
Illumina DMA Driver	4.5.3	4.5.3
Local Run Manager (LRM)	2.4.1	2.4.1

## I. MiniSeq System Suite v2.1.0

### NEW FEATURES:

- Minor updates to support new MiniSeq Control Software v2.1.0

## II. MiniSeq Control Software v2.1.0

### NEW FEATURES:

- Enable support of the new MiniSeq Rapid HO Reagent Kit (100 cys) on the MiniSeq System.
- Shortened to XY stage settle time.
- Optimization of paired-end read and cycle.

### IMPROVEMENTS:

- Resolved duplicated folder and file location issue from MiniSeq Control Software v2.0.